

The **V7P (and V7TP)** sensor transmitters contain a miniature **pressure sensor** which is recessed in the end cap. The V7P/TP was designed for *internal* implantation. Please discuss external attachment use on animals with VEMCO.

Please note the pressure sensor is covered with a **white** tacky silicone, therefore keep the V7P/TP free of debris before cleaning, and **do not** physically poke/push on the sensor. For more information, see your **Transmitter Specification Booklet**.



## Handling

- **Handle the V7P/TP tags very gently.** Use hands or rubber tipped forceps. Do not squeeze, bump or drop the tag. If you suspect that the tag has been subjected to harsh treatment, you should probably not use it to tag your fish. Record any relevant notes on handling.
- **Do not place the V7P/TP near another magnet.** This may adversely impact the function of the tag.

## Disinfecting

- **Do NOT use any caustic and acidic solutions or heat** to disinfect these transmitters. Once ready to disinfect the transmitter, please remove the yellow protective sticker covering the depth sensor.
- **VEMCO has tested both Betadine and Stanhexidine separately as disinfectants** on the V7P/TP transmitters, with no adverse effects to the depth sensor.



VEMCO Customer Support: (+1.902.450.1700)